

# **Blu10** Prestained Protein Ladder

03 JAN 2022

Catalog Number	Size	Range
PMB10-0500	500 μΙ	11 to 180 kDa

#### **Storage Conditions**

Stable for up to 2 weeks at 25°C.

Stable for up to 3 months at 4°C.

Stable for up to 24 months at -20°C.

## **Description**

The Blu10 Prestained Protein Ladder is a combination of 10 pre-stained proteins with molecular weights from 11 to 180 kDa. The 10 recombinant proteins are covalently coupled with blue chromophore, while 1 green band at 25 kDa and a red band at 75 kDa, serve as reference bands. The Blu10 Prestained Protein Ladder keeps track of the size and separation of proteins during SDS-polyacrylamide gel electrophoresis, approximating protein size, and validating Western transfer efficiency on PVDF, nylon, or nitrocellulose membranes.

#### Kit Content(s)

#### **Storage Buffer**

Approximately 0.2~0.4 mg/ml of each protein in buffer (20 mM Trisphosphate pH 7.5 at 25°C), 2% SDS, 0.2 mM Dithiothreitol, 3.6 M Urea, and 15% (v/v) Glycerol.

#### Required materials but not provided

- Vertical Electrophoresis system
- Power supplies
- Vortex or equivalent
- Microcentrifuge


## **Reaction Setup**

3  $\mu$ l or 5  $\mu$ l per loading for clear visualization during electrophoresis on 15-well or 10-well mini-gel, respectively. 2.5  $\mu$ l per well for general Western transferring.



# Guide for Molecular Weight Estimation (kDa)

Refer to the Blu10 Prestained Protein Ladder patterns in various electrophoresis conditions:

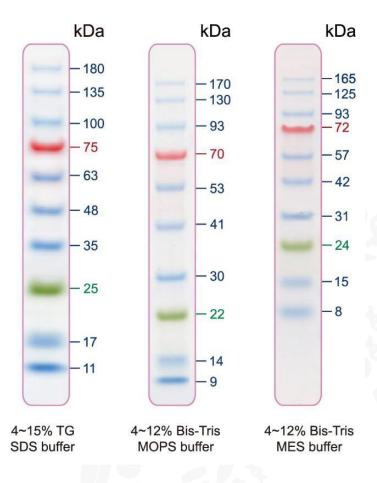

## Important notes

The molecular weight of each protein (kDa) was measured against an unstained protein ladder in every electrophoresis condition. Additional data should be considered for a more accurate adjustment.

